


PT3795-5

Cat. No. 632465

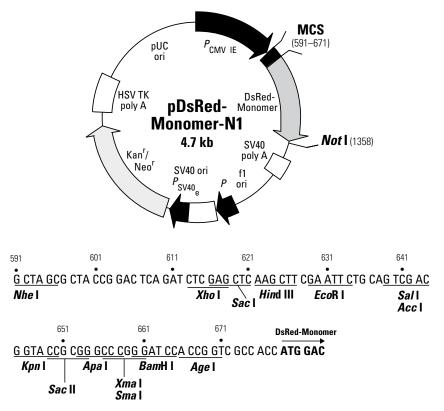

Restriction Map and Multiple Cloning Site (MCS) of pDsRed-Monomer-N1 Vector. Unique restriction sites are in bold. The *Not* I site follows the DsRed-Monomer stop codon. NOTE: The *Xba* I and *BcI* I sites are methylated in the DNA provided by Clontech Laboratories, Inc. If you wish to digest the vector with these enzymes, you will need to transform the vector into a dam– host and make fresh DNA.

Description

pDsRed-Monomer-N1 is a mammalian expression vector that encodes DsRed-Monomer (DsRed.M1), a monomeric mutant derived from the tetrameric *Discosoma* sp. red fluorescent protein DsRed (1). DsRed-Monomer contains forty-five amino acid substitutions (listed on page 2). When DsRed-Monomer is expressed in mammalian cell cultures, red fluorescent cells can be detected by either fluorescence microscopy or flow cytometry 12–16 hr after transfection (DsRed-Monomer excitation and emission maxima = 557 nm and 585 nm, respectively). The DsRed-Monomer coding sequence is human codon-optimized for high expression in mammalian cells (2).

DsRed-Monomer is well suited for use as a fusion tag. The multiple cloning site (MCS) in pDsRed-Monomer-N1 is positioned between the immediate early promoter of CMV ($P_{\text{CMV IE}}$) and the DsRed-Monomer coding sequence. Genes cloned into the MCS are expressed as fusions to the N-terminus of DsRed-Monomer if they are in the same reading frame as DsRed-Monomer and there are no intervening stop codons. A Kozak consensus sequence is located immediately upstream of the DsRed-Monomer gene to enhance translational efficiency in eukaryotic systems (3). SV40 polyadenylation signals downstream of the DsRed-Monomer gene direct proper processing of the 3' end of the DsRed-Monomer mRNA. The vector backbone contains an SV40 origin for replication in mammalian cells expressing the SV40T antigen, a pUC origin of replication for propagation in *E. coli*, and an f1 origin for single-stranded DNA production. A neomycin-resistance cassette (Neo¹) allows stably transfected eukaryotic cells to be selected using G418. This cassette consists of the SV40 early promoter, the neomycin/kanamycin resistance gene of Tn5, and polyadenylation signals from the Herpes simplex virus thymidine kinase (HSVTK) gene. A bacterial promoter upstream of the cassette confers kanamycin resistance in *E. coli*.

Use

pDsRed-Monomer-N1 can be used to construct fusions to the N-terminus of DsRed-Monomer. If a fusion construct retains the fluorescent properties of the native DsRed-Monomer protein, its expression can be monitored by flow cytometry and its localization *in vivo* can be determined by fluorescence microscopy. The target gene must be cloned into pDsRed-Monomer-N1 so that it is in frame with the DsRed-Monomer coding sequence, with no intervening in-frame stop codons. The inserted gene must include an initiating ATG codon. Recombinant pDsRed-Monomer-N1 can be transfected into mammalian cells using any standard transfection method. If required, stable transfectants can be selected using G418 (4). pDsRed-Monomer-N1 can also be used as a cotransfection marker; the unmodified vector will express DsRed-Monomer.

The DsRed1-N Sequencing Primer (Cat. No. 632387) can be used to sequence genes cloned adjacent to the 5' end of the DsRed-Monomer coding region.

For Western blotting, the Living Colors® DsRed Polyclonal Antibody (Cat. No. 632496) can be used to recognize the DsRed-Monomer protein. However, to generate optimal results it may be necessary to use a higher concentration of antibody than recommended on the DsRed Polyclonal Antibody Certificate of Analysis.

Location of features

Human cytomegalovirus (CMV) immediate early promoter: 1–589

Enhancer region: 59-465; TATA box: 554-560

Transcription start point: 583

C→G mutation to remove Sac I site: 569

• Multiple Cloning Site: 591-671

• Human codon-optimized DsRed-Monomer gene

Kozak consensus translation initiation site: 672–682 Start codon (ATG): 679–681; Stop codon: 1354–1356

Amino acid substitutions (DsRed→DsRed-Monomer)

GCC→GAC (Ala-2 to Asp) mutation: 682–684

TCC AAC (Ser-3 to Asn) mutation: 685–687

TCC→ACC (Ser-4 to Thr) mutation: 688–690

AAG→GAG (Lys-5 to Glu) mutation: 691–693

AAC→GAC (Asn-6 to Asp) mutation: 694–696

CGC→CAG (Arg-13 to Gln) mutation: 715–717

ACC→TCC (Thr-21 to Ser) mutation: 739–741

GAG→TAC (Glu-26 to Tyr) mutation: 754–756

CGC→AAG (Arg-36 to Lys) mutation: 784–786

CAC→ACC (His-41 to Thr) mutation: 799–801

AAC→CAG (Asn-42 to Gln) mutation: 802–804 GTG→GCC (Val-44 to Ala) mutation: 808–810

AAG -> CAG (Lys-47 to GIn) mutation: 817-819

GTG→GCC (Val-71to Ala) mutation: 889–891

AAG→ATG (Lys-83 to Met) mutation: 925–927

 $AAG \rightarrow ACC$ (Lys-92 to Thr) mutation: 952–954

 $GTG \rightarrow TCC$ (Val-96 to Ser) mutation: 964–966

ACC→GAG (Thr-106 to Glu) mutation: 994-996

ACC→CAG (Thr-108 to Gln) mutation: 1000–1002

TCC→ACC (Ser-117 to Thr) mutation: 1027–1029

ATC \rightarrow AAG (Ile-125 to Lys) mutation: 1051–1053

TCC→GCC (Ser-131 to Ala) mutation: 1069-1071

ATG→GCC (Met-141 to Ala) mutation: 1099-1101

GCC→CCC (Ala-145 to Pro) mutation: 1111–1113

CGC→AAG (Arg-149 to Lys) mutation: 1123-1125

CGC→CAG (Arg-153 to Gln) mutation: 1135-1137

CAC→TCC (His-162 to Ser) mutation: 1162–1164

AAG→CAC (Lys-163 to His) mutation: 1165–1167

CTG→ACC (Leu-174 to Thr) mutation: 1198–1200

GTG \rightarrow TGC (Val-175 to Cys) mutation: 1201–1203 GAG \rightarrow GAC (Glu-176 to Asp) mutation: 1204–1206

TCC→ACC (Ser-179 to Thr) mutation: 1213–1215

ATC \rightarrow GTG (Ile-180 to Val) mutation: 1216–1218

ATG-AAG (Met-182 to Lys) mutation: 1222-1224

TAC→AAC (Tyr-192 to Asn) mutation: 1252–1254

TAC \rightarrow CAC (Tyr-193 to His) mutation: 1255–1257 TCC \rightarrow AAC (Ser-203 to Asn) mutation: 1285–1287 ATC \rightarrow GTG (Ile-210 to Val) mutation: 1306–1308 CGC \rightarrow CAC (Arg-216 to His) mutation: 1324–1326 ACC \rightarrow GCC (Thr-217 to Ala) mutation: 1327–1329 GGC \rightarrow GCC (Gly-219 to Ala) mutation: 1333–1335 CAC \rightarrow TCC (His-222 to Ser) mutation: 1342–1344 CTG \rightarrow GGC (Leu-223 to Gly) mutation: 1345–1347 TTC \rightarrow TCC (Phe-224 to Ser) mutation: 1351–1353 CTG \rightarrow CAG (Leu-225 to Gln) mutation: 1351–1353

• SV40 early mRNA polyadenylation signal

Polyadenylation signals: 1510–1515 & 1539–1544; mRNA 3' ends: 1548 & 1560

- f1 single-strand DNA origin: 1607–2062 (Packages the noncoding strand of DsRed-Monomer)
- Bacterial promoter for expression of Kan^r gene:

-35 region: 2124-2129; -10 region: 2147-2152

Transcription start point: 2159

- SV40 origin of replication: 2403-2538
- SV40 early promoter

Enhancer (72-bp tandem repeats): 2236-2307 & 2308-2379

21-bp repeats: 2383-2403, 2404-2424 & 2426-2446

Early promoter element: 2459-2465

Major transcription start points: 2455, 2493, 2499 & 2504

Kanamycin/neomycin resistance gene

Neomycin phosphotransferase coding sequences: Start codon (ATG): 2587-2589; Stop codon: 3379-3381

G→A mutation to remove Pst I site: 2769

C→A (Arg to Ser) mutation to remove BssH II site: 3115

Herpes simplex virus (HSV) thymidine kinase (TK) polyadenylation signal

Polyadenylation signals: 3617–3622 & 3630–3635

pUC plasmid replication origin: 3966–4609

Sequencing primer location

DsRed1-N Sequencing Primer (Cat. No. 632387; 5'-GTACTGGAACTGGGGGGACAG-3'): 879–859

Propagation in E. coli

- Suitable host strains: DH5α, HB101 and other general purpose strains. Single-stranded DNA production requires
 a host containing an F plasmid, such as the JM109 or XL1-Blue strains.
- Selectable marker: plasmid confers resistance to kanamycin (50 µg/ml) in E. coli hosts.
- E. coli replication origin: pUC
- Copy number: high
- Plasmid incompatibility group: pMB1/ColE1

Excitation and emission maxima of DsRed-Monomer

- Excitation maximum = 557 nm
- Emission maximum = 585 nm

References

- 1. Matz, M. V., et al. (1999) Nature Biotech. 17:969-973.
- 2. Haas, J., et al. (1996) Curr. Biol. 6:315-324.
- 3. Kozak, M. (1987) Nucleic Acids Res. 15:8125-8148.
- 4. Gorman, C. (1985) In DNA cloning: A Practical Approach, Vol. II. Ed. D. M. Glover. (IRL Press, Oxford, U.K.), pp. 143-190.

Note: The attached sequence file has been compiled from information in the sequence databases, published literature, and other sources, together with partial sequences obtained by Clontech Laboratories, Inc.. This vector has not been completely sequenced.

Notice to Purchaser

Clontech products are to be used for research purposes only. They may not be used for any other purpose, including, but not limited to, use in drugs, *in vitro* diagnostic purposes, therapeutics, or in humans. Clontech products may not be transferred to third parties, resold, modified for resale, or used to manufacture commercial products or to provide a service to third parties without written approval of Clontech Laboratories, Inc.

Not-For-Profit Entities: Orders may be placed in the normal manner by contacting your local representative or Clontech Customer Service at 650.919.7300. At its discretion, Clontech grants Not-For-Profit Entities a non-exclusive, personal, limited license to use this product for non-commercial life science research use only. Such license specifically excludes the right to sell or otherwise transfer this product, its components or derivatives thereof to third parties. No modifications to the protein coding sequence may be made without express written permission from Clontech. Any other use of this product requires a license from Clontech. For license information, please contact a licensing representative by phone at 650.919.7320 or by e-mail at licensing@clontech.com.

For-Profit Entities wishing to use this product are required to obtain a license from Clontech. For license information, please contact a licensing representative by phone at 650.919.7320 or by e-mail at licensing@clontech.com.

The CMV promoter is covered under U.S. Patent Nos. 5,168,062, and 5,385,839 assigned to the University of Iowa Research Foundation. This product is covered under U.S. Patent No. 7,250,298.

Clontech, the Clontech logo and all other trademarks are the property of Clontech Laboratories, Inc., unless noted otherwise. Clontech is a Takara Bio Company. ©2008 Clontech Laboratories, Inc.

Clontech Laboratories, Inc.

www.clontech.com

Protocol No. PT3795-5

Version No. PR842501