

Restriction map of pRevTet-Off<sup>™</sup> Vector. Unique restriction sites are in bold.

### **Description:**

pRevTet-Off is a retroviral vector expressing the tetracycline-controlled transactivator (tTA) from the CMV promoter. This vector is derived from pLNCX, a retroviral vector created using elements of Moloney murine leukemia virus (MoMuLV) and Moloney murine sarcoma virus (MoMuSV) as described (1). tTA is a fusion of amino acids 1–207 of the tet repressor (TetR) and the negatively charged C-terminal activation domain (130 amino acids) of the VP16 protein of Herpes Simplex Virus. The 5' viral LTR controls expression of the transcript that contains  $\Psi^+$  (the extended viral packaging signal), and the neomycin resistance gene (Neo<sup>r</sup>) for antibiotic selection in mammalian cells. tTA is derived from vectors described previously (2, 3). pRevTet-Off also includes the *E. coli* Amp<sup>r</sup> gene for antibiotic selection in bacteria.

### Use:

pRevTet-Off can be used to establish stable Tet-Off cell lines via retrovirus-mediated gene transfer (4). Retroviral gene transfer allows the highly efficient transduction of virtually all dividing cell types. The RevTet Systems are also suitable for establishing transgenic animals. In combination with the pRev-TRE retroviral expression vector, a gene of interest can be inducibly expressed at high levels in response to varying concentrations of tetracycline (Tc) or Tc derivatives such as doxycycline (Dox). tTA binds to the Tet-response element (TRE), thus activating transcription in the absence of Tc or Dox. AsTc or Dox is added to the culture medium, transcription from the inducible promoter is turned off in a highly dose-dependent manner. pRevTet-Off lacks the viral genes *gag*, *pol*, and *env*, which are supplied by the packaging cell line. It can be transfected into a high titer packaging cell line and thereby mediate production of infectious, replication-incompetent retroviral particles (1, 5–6). The transcript produced by the pRevTet-Off construct is recognized by the viral structural proteins expressed in a packaging cell line and packaged into infectious retroviral particles. Because the RNA transcript packaged in these particles does not contain the viral genes, it cannot replicate in the target cells that it infects.

The level of induction in cell populations infected with this vector depends on the efficiency of infection, the site of integration, and the titer of the virus. Viral supernatants with titers  $>10^5$  cfu/ml should be produced to achieve high-level induction.

## Location of features:

- 5' MoMuSV LTR: 1-589
- $\Psi^+$  (extended packaging signal): 659–1468
- Neomycin resistance gene: 1512–2306
  - Start codon: 1512–1514; stop codon: 2304–2306
- CMV promoter (P<sub>CMV</sub>): 2655–3465
- tTA coding sequence: 3476-4483
- 3' MoMuLV LTR: 4749-5342
- Ampicillin resistance (β-lactamase) gene: 7498–6638
  Start codon: 7498–7495; stop codon: 6640–6638

# Propagation in *E. coli*:

- Suitable host strains: DH5α, HB101, and other general purpose strains.
- Selectable marker: plasmid confers resistance to ampicillin (50 µg/ml) to E. coli hosts.
- E. coli replication origin: Col E1
- Copy number: low copy

### **References:**

- 1. Miller, A. D. & Rosman, G. J. (1989) *BioTechniques* **7**:980–990.
- 2. Tet Expression Systems and Cell Lines (July 1996) *Clontechniques* XI(3):2–5.
- 3. Gossen, M. & Bujard, H. (1992) Proc. Natl. Acad. Sci. USA 89:5547-5551.
- 4. RevTet-On & RevTet-Off Systems (April 1998) Clontechniques XIII(2):8-9.
- 5. *Retroviruses* (1996) CSHL Press.
- 6. Current Protocols in Molecular Biology (1997) Supplement 36, Section III

### Notes:

The attached sequence file has been compiled from information in the sequence databases, published literature, and other sources, together with partial sequences obtained by Clontech. This vector has not been completely sequenced.

The viral supernatants produced by this retroviral vector could, depending on your cloned insert, contain potentially hazardous recombinant virus. Due caution must be exercised in the production and handling of recombinant retrovirus. Appropriate NIH, regional, and institutional guidelines apply.

### Notice to Purchaser

Clontech products are to be used for research purposes only. They may not be used for any other purpose, including, but not limited to, use in drugs, *in vitro* diagnostic purposes, therapeutics, or in humans. Clontech products may not be transferred to third parties, resold, modified for resale, or used to manufacture commercial products or to provide a service to third parties without written approval of Clontech Laboratories, Inc.

Retroviral Vectors are sold under license from the Fred Hutchinson Cancer Research Center. Rights to use this product are limited to research only. No other rights are conveyed. Inquiry into the availability of a license to broader rights or the use of this product for commercial purposes should be directed to Fred Hutchinson Cancer Research Center, Technology Transfer Office, 1100 Fairview Avenue North, J6-200, Seattle, WA 98109. Purchase of this product does not grant rights to: (1) offer the materials or any derivatives thereof for resale; or (2) to distribute or transfer the materials or any derivatives thereof to third parties.

Use of the Tetracycline controllable expression systems (the "Tet Technology") is covered by a series of patents including U.S. Patent Nos. 5,464,758 and 5,814,618, which are proprietary toTET Systems Holding GmbH & Co. KG. Academic research institutions are granted an automatic license with the purchase of this product to use the Tet Technology only for internal, academic research purposes, which license specifically excludes the right to sell, or otherwise transfer, the Tet Technology or its component parts to third parties. Notwith-standing the above, academic and not-for profit research institutions whose research using the Tet Technology is sponsored by for profit organizations, which shall receive ownership to all data and results stemming from the sponsored research, shall need a commercial license agreement from IP Merchandisers in order to use the Tet Technology. In accepting this license, all users acknowledge that the Tet Technology is experimental in nature. TET Systems Holding GmbH & Co. KG makes no warranties, express or implied or of any kind, and hereby disclaims any warranties, representations, or guarantees of any kind as to the TetTechnology, patents, or products. All others are invited to request a license from TET Systems Holding GmbH & Co. KG prior to purchasing these reagents or using them for any purpose. Clontech is required by its licensing agreement to submit a report of all purchasers of the Tet-controllable expression system

Hans Peter Kneubuehl TET Systems Holding GmbH & Co. KG Im Neuenheimer Feld 582 69120 Heidelberg Germany Tel +49 6221 588 04 00 Fax +49 6221 588 04 04 eMail: kneubuehl@tet-systems.de or use our electronic licensing request form via http://www.tetsystems.com/main\_inguiry.htm

Clontech, the Clontech logo and all other trademarks are the property of Clontech Laboratories, Inc. Clontech is a Takara Bio Company. ©2006